DOC022.98.00648

TOC-X5

10/2021, Edition 4

Ръководство за потребителя	bg
Základní uživatelská příručka	cs
Grundlæggende brugervejledning	da
Bedienungsanleitung	de
Βασικό Εγχειρίδιο Χρήστη	el
User Manual	en
Manual del usuario	es
Basic käyttöohjeet	fi
Manuel d'utilisation	fr
Osnovni korisnički priručnik	hr
Alap felhasználói kézikönyv	hu
Manuale utente di base	it
Basisgebruikershandleiding	nl
Podstawowa instrukcja obsługi	pl
Manual do utilizador básico	pt
Manual al utilizatorului	ro
Основное руководство пользователя	ru
Základný návod na použitie	sk
Osnovni uporabniški priročnik	sl
Grundläggande bruksanvisning	sv
Temel Kullanıcı Kılavuzu	tr

8Български	3
Čeština	9
Dansk	15
Deutsch	21
Ελληνικά	27
English	33
Español	39
Suomi	45
Français	51
Hrvatski	57
Magyar	63
Italiano	69
Nederlands	75
Polski	81
Português	87
Română	93
Русский	99
Slovenský jazyk	105
Slovenščina	111
Svenska	117
Türkçe	123

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Specifications

Specifications are subject to change without notice.

Specification	Details
Dimensions	190 x 112 x 240 mm (7,48 x 4,41 x 9,45 in.)
Power supply	100–240 VAC $\pm$ 10%; 50/60 Hz; 0,75 A
Instrument	15 VDC; 30 VA
Plug-in power supply unit	LZV563
Operating requirements	10 to 40 °C (50 to 104 °F); maximum 80% relative humidity (non condensing)
Storage requirements	-40 to 60 °C (-40 to 140 °F); maximum 80% relative humidity (non condensing)
Weight	2,9 kg (6,39 lb), inclusive plug-in power supply unit
Enclosure	IP20
Overvoltage category	II
Pollution degree	2
Altitude	2000 m (6562 ft) maximum
Environmental conditions	Indoor use
Compliance	CE
Warranty	EU: 2 years (US: 1 year)

General information

In no event will the manufacturer be liable for direct, indirect, special, incidental or consequential damages resulting from any defect or omission in this manual. The manufacturer reserves the right to make changes in this manual and the products it describes at any time, without notice or obligation. Revised editions are found on the manufacturer's website.

Safety information

The manufacturer is not responsible for any damages due to misapplication or misuse of this product including, without limitation, direct, incidental and consequential damages, and disclaims such damages to the full extent permitted under applicable law. The user is solely responsible to identify critical application risks and install appropriate mechanisms to protect processes during a possible equipment malfunction.

Please read this entire manual before unpacking, setting up or operating this equipment. Pay attention to all danger and caution statements. Failure to do so could result in serious injury to the operator or damage to the equipment.

Make sure that the protection provided by this equipment is not impaired, do not use or install this equipment in any manner other than that specified in this manual.

Use of hazard information

 **DANGER**

Indicates a potentially or imminently hazardous situation which, if not avoided, will result in death or serious injury.

 **WARNING**

Indicates a potentially or imminently hazardous situation which, if not avoided, could result in death or serious injury.

 **CAUTION**

Indicates a potentially hazardous situation that may result in minor or moderate injury.

NOTICE

Indicates a situation which, if not avoided, may cause damage to the instrument. Information that requires special emphasis.

Precautionary labels


Read all labels and tags attached to the instrument. Personal injury or damage to the instrument could occur if not observed. A symbol on the instrument is referenced in the manual with a precautionary statement.

	This symbol, if noted on the instrument, references the instructions manual for operation and/or safety information.
	Electrical equipment marked with this symbol may not be disposed of in European public disposal systems. Return old or end-of-life equipment to the manufacturer for disposal at no charge to the user.

Product overview

The TOC-X5 shaker is used in laboratories to expel total inorganic carbon (with the TOC cuvette tests LCK 385 / 386 / 387). The vibration unit can shake up to 8 vials simultaneously. The ports are labeled 1-8. The cycle time of a shaking process is 5 minutes. Refer to [Figure 1](#).

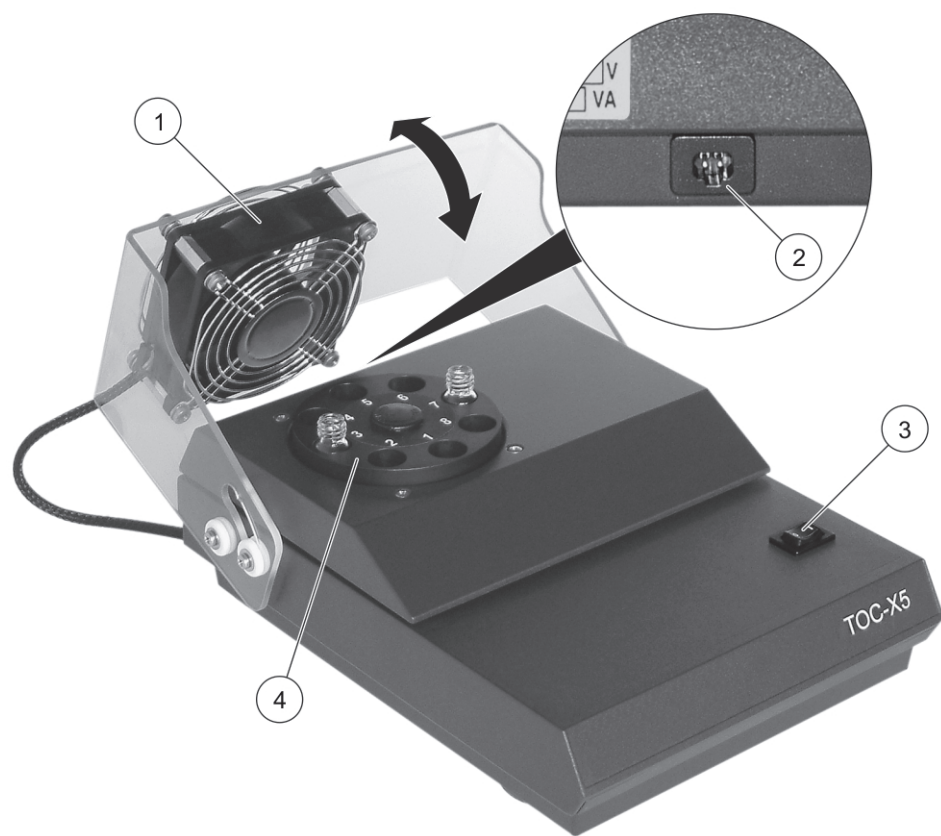Figure 1 Product overview

1	Pivoted safety cover with fan	3	Power switch
2	Interface for plug-in power supply unit	4	Vibration unit

Product components

Make sure that all components have been received. If any items are missing or damaged, contact the manufacturer or a sales representative immediately.

Shipped items:

- TOC-X5 shaker
- Plug-in power supply unit
- Non-slip mat

Installation guideline

DANGER

Multiple hazards. Only qualified personnel must conduct the tasks described in this section of the document.

Install the instrument:

- In a clean, dry, level, well ventilated, temperature controlled location
- In a location where the power switch and power socket are visible and easily accessible because the plug-in power supply is used as a separator in the case of a hazard

The instrument does not move with the suction cups installed. Make sure that the suction cups are clean and free from damage. Use the supplied non-slip mat, if the suction cups lose their adhesion on rough surfaces.

Operation

WARNING

Pivot the safety cover with fan over the vibration unit before the instrument is in use.

WARNING

**Do not insert fingers or any other objects through the protective grid of the fan.
Take care: Long hair can be sucked into the fan by the air current.**

NOTICE

Use only the power supply that is specified for this instrument.

NOTICE

**Do not grasp on or in the vibration unit during the shaking procedure.
Do not install a vial during the shaking procedure.
Do not remove a vial during the shaking procedure.
Set the power switch to off and wait until the vibration unit and the fan to stop before any activity.**

Obey these safety precautions:

- Do not use damaged glassware or apparatus. Discard all damaged equipment and replace it.
- Make sure that the vials are filled up to maximal 3/4

Always do the steps that follow in the order shown:

1. Connect the plug-in power supply unit to a power socket.
2. Install the vials fully into the vibration unit (up to ground contact).
3. Pivot the safety cover with fan over the vibration unit.
4. Set the power switch to on.
The shaking procedure starts. The timer will beep after five minutes and the procedure stops.
5. Pivot the safety cover with fan to the back.
6. Remove the vials.
7. Disconnect the plug-in power supply unit from the power socket.

The instrument can be used for shorter shaking times. Set the power switch to off any time to interrupt the shaking procedure.

Maintenance

DANGER

Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.

CAUTION

Multiple hazards. Only the manufacturer or an authorized representative must conduct all service activities. Do not disassemble the instrument for maintenance. If the internal components must be cleaned or repaired, contact the manufacturer.

CAUTION

Personal injury hazard. Broken glass can cause cuts. Use tools and personal protective equipment to remove broken glass.

Clean the instrument

Keep the instrument clean to get continuous and accurate operation.

NOTICE

Never use cleaning agents such as turpentine, acetone or similar products to clean the instrument.

NOTICE

Do not leave visible moisture on the instrument. Moisture can penetrate and damage the electronics inside.

1. Set the instrument off and disconnect the plug-in power supply unit.
2. Clean the surface of the instrument with a soft, moist cloth and a weak soap solution.
3. Dry the surface of the instrument with a lint-free cloth.

How to clean spills

CAUTION

Chemical exposure hazard. Dispose of chemicals and wastes in accordance with local, regional and national regulations.

1. Obey all facility safety protocols for spill control.
2. Discard the waste according to applicable regulations.

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P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
(800) 227-4224 (U.S.A. only)